

CertoClav Classic

The economical end-to-end solution made of stainless steel

The robust CertoClav Classic is the result of the further development of the EL made of stainless steel and is extremely low maintenance. A precision thermometer, a control valve, and a precision manometer with drag indicator provide precise processes and accurate results for all applications. This model is available in two versions, each with two temperature levels.

You can choose between 125/140 °C and 115/121 °C. These temperature settings are controlled via the control valve on the lid. These valves are not interchangeable, because the thermostat of the device is harmonized and calibrated with the respective valve!

A timer is integrated to regulate the duration of the sterilization process. If the timer is activated, the device switches off automatically after the sterilization time has passed. If the timer has not been activated, the autoclave must be switched off manually.



control valve

Important features

- robust construction
- 2 temperature levels
- simple operation
- integrated timer function
- venting by flow procedure
- precision thermometer
- precision manometer with drag pointer
- heating plate integrated into base
- thermostat controlled

Scope of delivery:

incl. instrument tray and power cable

Specifications

	CertoClav Classic 125/140 °C	CertoClav Classic 115/121 °C
Article number	8510174	8510184
Chamber volume	18 l	18 l
Chamber diameter	26 cm	26 cm
Chamber depth	35 cm	35 cm
Chamber material	stainless steel 316L	stainless steel 316L
Max. operating pressure	2,7 bar	2,7 bar
Mains connection	230 V, 50/60 Hz	230 V, 50/60 Hz
Heating capacity	1300 W	1300 W
Max. sterilization temperature	125 °C oder 140 °C	115 °C oder 121 °C
Net weight	8,8 kg	8,8 kg
Product dimensions (L*W*H)	44,5*30*60,5 cm	44,5*30*60,5 cm

Accessories

Article	Article number
Wire basket with handle	1300149
Service Set 2	1310186
Exhaust steam container	8582020
Exhaust steam hose	8582016
Distiller for water	8500482