

Pipette washer PSD



PSD is an easy to use electronic pipette washer dedicated to clean all the glass pipettes generally used.

It features a fixed program cycle including automatic addition of detergent which guarantees constant quality of the cleaning results. Soaking the pipettes in the heated solution of *VE-water and detergent will cause all the dirt particles to diffuse into the cleaning solution. The following rinsing cycles with tap water take away all the dirt particles out of the tank. Afterwards rinsing cycles with *VE-water avoid adhesion of lime of the pipettes thoroughly. Heating up the empty cleaning tank slowly will dry the pipettes in a smooth way in order to have them ready to use at the end of the program.



Technical information

Connection

Voltage: 230V / 16 A Current: 3,0 KW

Tap water: 3/4"; max. 60°C intake temperature

*VE-water: 3/4"

Water pressure: between 0,5 and 6 bar

Waste water: **DN20**

Consumption

lowhigh water level 4.9 kWh 7 kWh Energy:

*VE-water: max. 90 liter max. 130 liter Tap water: max. 180 liter max. 270 liter 0-400 ml 0-400 ml Detergent:

Temperature range

Cleaning: 30°C or 96°C 130°C Drying:

Programs for two kinds of water

Number: 1

Duration: ca. 8 hours

Capacity: max. 9 baskets

Weight/ material: 50 kg / stainless steel housing

Attachement: 8 kg

Dimensions:

breadth: 325 mm depth: 530 mm

height: 920 mm (device top edge with closed cover plate)

height: 1230 mm (with open cover plate)

Cleaning capacity

| Messure pipettes | | | | | | | |
|------------------|--------|--------|--------|---------|---------|---------|----------|
| Volume | 1,0 ml | 2,0 ml | 5,0 ml | 10,0 ml | 20,0 ml | 50,0 ml | 100,0 ml |
| Per rinser | 1305 | 1080 | 810 | 441 | 324 | 144 | 54 |
| Full pipettes | | | | | | | |
| Volume | 2,0 ml | 3,0 ml | 5,0 ml | 10,0 ml | 20,0 ml | 50,0 ml | 100,0 ml |
| Per rinser | 900 | 576 | 324 | 225 | 144 | 54 | 20 |



Equipment

6 pcs. Baskets 80x80x415 mm 3 pcs. baskets 80x80x300 mm 9 pcs. Numbering clambs

to mark the baskets

5 liter liquid purifying agent

Technical documentation and

connection set

1 pcs. connection set:

2x water inlet house 1x water outlet house

1x hose bow 90° 6x hose clamp 4x hose coupling

4x seal

Stand: 12.04.2012